Solexa Data Crunching

David Lapointe March 20, 2008


Data Crunching

- Binar A Tour
- Resources on Binar
- How Data is Crunched
- Other Resources

Binar



Binar: Components

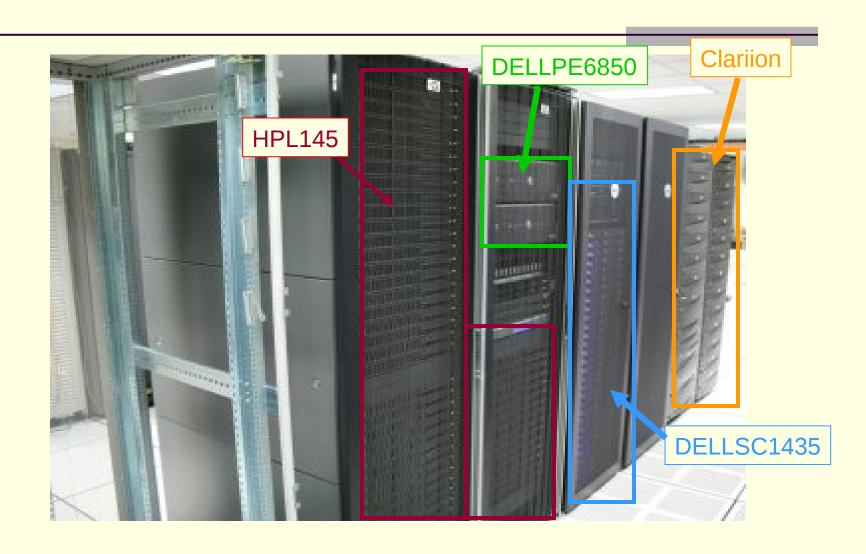

More Specs

- HPL145
 - 2 dual core Opteron/4 Gb mem
 - 55 nodes
- DELLPE6850
 - 4 quad core Intel/64 Gb mem
 - 2 nodes
- DELLSC1435
 - 2 dual core Opteron/4 Gb mem
 - 25 nodes
 - Infiniband capable

Cluster Use/Etiquette

- Use queue system to run jobs
 - Sun Grid Engine queueing
 - Fair share queueing is in use
- File space is not storage
 - Do not store data on the cluster
 - Cluster has 2.0 Tb only for workspace
 - 23 Active Users

Queues

\$qsub script basic form

\$qstat shows jobs in queue

\$qstat | grep user shows jobs for username

Info for Binar and Sun Grid Engine is online at http://inside.umassmed.edu/is/acs/ResearchComputing/researchclusters.aspx

Script Template for Queue

```
#!/bin/bash
#$ -S /bin/bash <- specify shell
                   <- use current working dir
#$ -cwd
#$ ... other sge parameters
PATH=<path to your program>:$PATH
export PATH
...rest of script ....
Then submit script within working directory
$qsub -cwd script
```

Scripts

It is very important to set custom paths, esp for programs that you write, or using bioperl.

/share/... is common to all nodes.
/export/home/<user> (your home directory) is available to remote nodes.

Important Directories

/share/apps/bin commonly used programs /share/apps/pipeline/... Solexa pipeline progs

/share/nemo/Genomes contains genome directories formatted for Eland

/share/nemo/Genomes/hg18

/share/nemo/Genomes/mm9

Resources on Binar

- EMBOSS (similar to GCG, scriptable)
- Clustalw, T_Coffee multiple alignment
- Phylip, MrBayes phylogenetic
- BLAST, mpi-blast
- HMMER
- Bioperl, biopython

These have their own directories.

Current Genomes on Binar

hg18 Human (UCSC)

mm9, mm8 Mouse (UCSC)

dm5.5, dm5.4 Drosophila (Flybase)

yeast SGD

z7v
Zebrafish

ceWS187 C.Elegans

/share/nemo/Genomes/xxxx

Creating a Genome file for Eland

You can run eland against a custom Genome.

- 2. Each piece (e.g. a chromosome) must be a separate fasta file.
- Create a directory for the genome mkdir ~/mygenome
- 4. Run squashGenome /share/nemo/pileline/Eland/squashGenome \ ~/mygenome pathtoFasta/*.fasta
- 5. More detail in Pipeline docs

Pipeline processing

- Three phases
 - Image Analysis -Firecrest
 - Base Calling Bustard
 - Sequence Mapping —Gerald
- Initial Run to generate Cross-talk and offsets
- Run the first two by lane
 - If you specify Genome Gerald can be run
 - If we don't have the genome, let us know what and where, then we can install it.

Post-processing

- When runs are done several files are available
 - quality files, base calls
 - sequences, remapping results
 - summary data
- These are packaged and delivered to you.
- Let me know if you need custom programs for analysis, i.e. ones not delivered with the pipeline

Resources

- Cluster info http://inside.umassmed.edu/is/acs/ResearchComputing/researchclusters.aspx
- Pipeline Info http://biotools.umassmed.edu/BIOCORE/pipeline
- Solexa Google Group http://groups.google.com/group/solexa
- Bioc-Seq (New bioconductor group) https://stat.ethz.ch/mailman/listinfo/bioc-sig-sequencing
- UNIX On-line Help http://biotools.umassmed.edu/unixhelp